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CATALOG

Laboratory equipment for environmental
control of water, soil and air

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Sampling Equipment
Chromatography
Organic Compounds Analysers
Equipment for air quality monitoring
Measuring equipment for meteorological
parameters

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Dear partners!

Topan LLP is a company, which provides services and equipment delivery for industrial enterprises of Kazakhstan.

One of the direction of activities of Topan LLP is the organization and complex equipping of environmental monitoring posts for oil and gas companies (deposits, oil and gas refineries), for the chemical industry, steel plants, cement plants, power plants and water treatment facilities (with optional equipment for water analysis).

In our catalog we offer you to explore the wide range of **equipment for ecological control of water, air and soil** from the leading manufacturers of these products.

For more detailed information please contact us or send your request to our email address.

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CONTENT:

Sampling Equipment	
Liquid Sampling Devices	4
Soil Sampling Devices	5
General Physico-Chemical Parameters	
Turbidity and Color Analyzers	6
Laboratory Meters; pH, ISE, conductivity	10
Automatic Titrators	11
Vacuum Filtration Apparatuses	12
Mass measurements (analytical balances)	13
Ion Chromatography	15
Atomic Absorption Spectroscopy	16
UV-Vis-Spectroscopy	17
High Performance Liquid Chromatography	21
Organic Compounds Analysers	22
Analysers of Oil products in Water and Soils	23
Total Organic Carbon (TOC) of Water and Soil	24
Equipment for air quality monitoring	
Stationary gas analysers	25
Equipment for calibration of gas analysers	28
Portable gas analysers	30
High Volume Samplers	33
Measuring equipment for meteorological parameters	
Wind speed and Wind direction sensors	34
Ambient temperature, relative humidity and barometric pressure sensors	35
Brief operation description of AQMS (air quality monitoring station)	37

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LIQUID SAMPLING DEVICES

Dip Sampler, plunging siphon (Stechheber)

Dip sampler for simple removal of liquid samples. Good view due to transparent tubes. Upper end with reflux valve, lower end with ball valve.

- PP, transparent
- Dip sampler, steplessly extendable from 100 cm to 200 cm
- All-layer and point sampling possible

Use

All-layer sample

1. Insert the dip sampler slowly into the liquid with the valve open.
2. Close valve.
3. Pull the dip sampler back and take out the sample.

Point sampling/soil sample

1. Insert the dip sampler to the targeted point/to the ground with closed valve.
2. Open the valve - the liquid flows in.
3. Close the valve.
4. Pull out the dip sampler, take out the sample.

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ChemoSampler

The ChemoSampler was developed for sampling of aggressive liquids, e.g. acids, alkalis, detergents. The samples are drawn directly into the screw-on original sampling bottle. The tamperproof bottle closure of these bottles with original safety ring ensures unadulterated sample transport. Samples can be taken from depths of as much as 200 cm, e.g. from tanks, silos or tankers, by using the extension rod.

- PP, grey
- Resistant to acids and alkalis
- Incl. sample bottle 1000 ml
- The set includes 1000 ml sample bottle and 100 cm extension rod

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SOIL SAMPLING DEVICES

Mole with handle and bar

Very robust model with specially hardened forged drill bits. For determining soil properties, moisture and structure. The shape of the Mole's drill bits ensures minimum friction when drilling and removing the Mole from the ground. This reduces physical effort to a minimum.

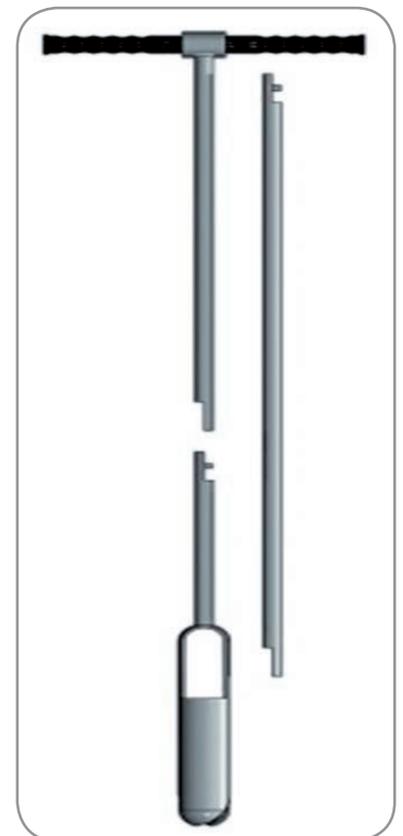
The Mole has different components:

- handle with 75 cm rod;
- extension rod 100 cm if necessary;
- and one of seven available drill bits, each of which is designed for a particular soil type.

There are various drill bits for Mole, which are each designed for very specific soil properties:

- Clay and damp soil;
- Sand;
- Coarse sand;
- Combination;
- Riverside;
- Gravel drill;
- Stonecatcher.

The components can be combined without any problems. Extension rods can extend the working depth of the Mole down to 5 m.

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TURBIDITY AND COLOR ANALYZERS

Turbidity can be interpreted as a measure of the relative clarity of water. Turbidity is not a direct measure of suspended particles in water but, instead, a measure of the scattering effect such particles have on light.

Today, many methods exist for the determination of water contaminants, yet turbidity measurement is still important because it is a simple and undeniable indicator of water quality change. A sudden change in turbidity may indicate an additional pollution source (biological, organic or inorganic) or may signal a problem in the water treatment process.

2100 series turbidimeters are designed for laboratory measurements of turbidity with high accuracy, and for the express control on the production line.

- laboratory turbidimeters 2100N, 2100N IS, 2100AN, 2100AN IS
- portable turbidimeters 2100Q/2100Q IS

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2100 Series Laboratory Turbidimeters (2100N, 2100N IS, 2100AN, 2100AN IS)

2100 Series Laboratory Turbidimeters are engineered to provide superior accuracy and sensitivity in any application. The 2100N Laboratory Turbidimeter is equipped with a stable halogen-filled tungsten filament lamp.

- **2100N/2100N IS** - **N** Series for standard applications.
- **2100AN/2100AN IS** - **AN** Series for professional applications.

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Portable Turbidimeters 2100Q/2100Q IS

2100Q and 2100Q IS - portable turbidimeter offer unrivaled ease of use and accuracy of measurement of turbidity. They can be used both in laboratory and in field studies.

- Easy on-screen assisted calibration and verification
- Simple data transfer
- Accurate for rapidly settling samples
- Convenient data logging
- Optical system for precision in the field

2100Q Turbidimeter meets the requirements of USEPA Method 180.1 with an incandescent lamp.

2100Q IS Turbidimeter meets the requirements of ISO 7027 with an infrared LED.

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Specification	2100Q	2100Q IS
Measurement Modes	NTU	FNU
Regulatory	EPA 180.1	ISO 7027
Accuracy	±2% (0 – 1000 NTU)	
Repeatability	±1%; ±0.01% NTU	
Resolution	0.01	0.01
Light Source	incandescent lamp.	infrared LED
Source	4 AA batteries	

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Color of Water Portable Photometer HI 96727.

Basic features:

- measures the true and apparent color in water and wastewater in the 0 to 500 PCU;
- uses an exclusive positive-locking system to ensure that the cuvette is in the same position every time it is placed into the measurement cell;
- designed to fit a cuvette with a larger neck making it easier to add both sample and reagents;
- the cuvette is made from special optical glass to obtain best results.

Operating conditions:

- Ambient temperature 0 ... 50 ° C;
- Relative humidity up to 95% (without condensation).

Standard delivery set:

- HI 96727 Colorimeter with instruction manual
- sample cuvettes with caps (2) CAL CHECK™
- Certified calibration and verification standards
- 9V battery and BEPS (Battery Error Prevention System)

Specification

- Range: 0 to 500 PCU (Platinum Cobalt Units)
- Resolution: 10 PCU
- Accuracy: ± 10 PCU $\pm 5\%$ of reading @ 25°С
- Light Source: tungsten lamp
- Light Detection: silicon photocell with narrow band interference filter @ 420 nm
- Battery Type: Life 9V approx. 40 hours of continuous use; auto-off after 10 minutes of non use

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LABORATORY METERS; PH, ISE, CONDUCTIVITY

Seven Compact Series and Seven Excellence Series Laboratory Multi-parameter Meters

The **S220 SevenCompact™** pH/Ion universal well designed meter is the perfect choice for a wide range of applications. It is much more than a pH meter: it can also measure ORP and ion concentration in various units.

The included electrode holder with the perfectly vertical uPlace™ movement helps you bring your electrode in the position where it works best for your sample. This allows for faster measurements and reduces the risk of the sample vessel tipping over and/or damage the sensor.

SevenCompact™ gives you the choice to view all measurement information on the display or to focus on the essentials with the uFocus™ view. Switching between the different display layouts is easy, requiring only a single keypress.

The SevenExcellence multi-parameter bench meter for precise measurements of pH, conductivity, ion concentration and ORP (Oxidation Reduction Potential).

Measuring range:

-2.000 ... 20.000 pH with selectable discreteness: 0.001 / 0.01 / 0.1;

-1999.9 ... 1999.9 mV;

0 ... 999,999 g/l;

-2 ... 20 pX;

temperature -30 ... 130 C;

manual or automatic temperature compensation;

3 options of recording the results of measurements: automatic, manual or time based.

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АНИОН Series Laboratory Single-parameter and Multi-parameter Meters

- Simplicity and ease of use;
- High technical and technological level;
- -All necessary metrological characteristics;
- -Developed set of service functions
- High quality of manufacture;
- -Affordable prices.

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Titration Excellence series titrators are recommended for analysis by titration with potentiometric, voltammetric and photometric indication. Types of analysis: chelatometry, turbidimetry, acid-base titration, the two-phase titration of surfactants and other

For the purposes of determining the total water hardness, titrators are equipped with a sensor for color transition titration and turbidimetric titration – DP5 Fototrod.

Five wavelengths (manual switch): 520, 555, 590, 620 and 660 nm.

One Click Titration™ – Intuitive and easy to understand

- Unsurpassed user guidance
- Secure navigation
- Easily selectable and adaptable methods

Improved efficiency – Through intelligence and automation

- One Click Titration™
- Series sequence
- Task button

Complete overview and security – Thanks to comprehensive control

- Intelligent burettes
- Automatic recognition of burette drives and sensor boards
- Dual Mode

For standard applications

Everything you need for titration, delivered as one kit. Besides the main module, touch terminal and burettes a kit may include automatic feeder and the titration stand.

You can also order:

- electrodes and sensors, changeable burettes and accessories (depending on the titration method, must be ordered separately);
- the ability to upgrade by three dosing modules (for addition of reagents), to replace the pH board on conductivity board;
- Professional management of user rights;
- Interfaces: USB, Ethernet, RS232, CAN.

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Vacuum Filtration Apparatuses

Purpose:

- Vacuum filtration of drinking water samples in sanitary and microbiological analyzes in accordance with GOST 18963-73, SanPin 2.1.4.1074-01, SanPin 2.1.4.1116-02 and MUK 4.2.1018-01;
- Microbiological testing in food, medical, pharmaceutical and other industries;
- Fine filtration of water in physical and chemical laboratories and research projects;
- Analysis of water to determine the content of suspended solids and dry residue of water;
- Determination of the purity of petroleum products, fuels and oils;
- Turbidity and water color;
- Preparation of samples in physico-chemical analysis of water.

To perform specific tasks there are different modifications of devices:

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MASS MEASUREMENTS (ANALYTICAL BALANCES)

Sartogosc CE Series Analytical Balances

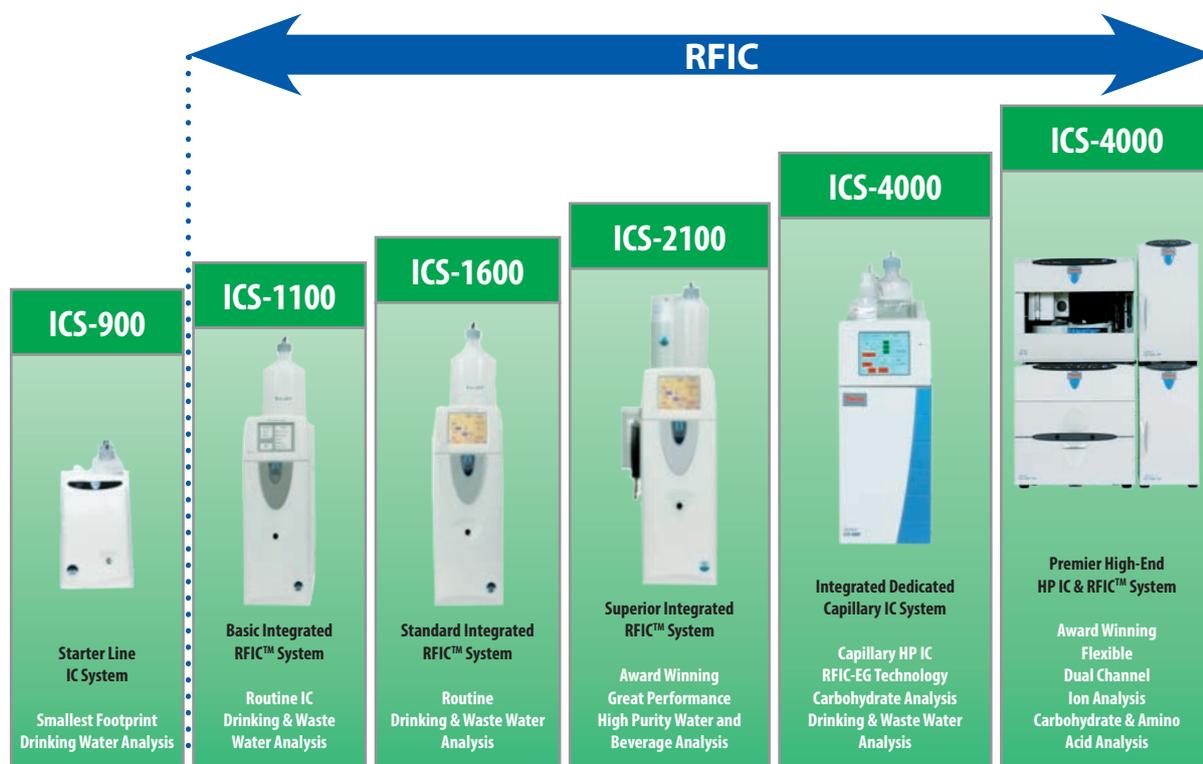
CE Series analytical balances are reliable and easy to operate, equipped with a wide range of applications and are indispensable in:

- laboratories;
- controlling organizations;
- pharmaceuticals;
- educational and research projects;
- in the production process.

Benefits:

- Expanded product range;
- Conform to international quality standards;
- Selection of variants;
- Text Menu in Russian.

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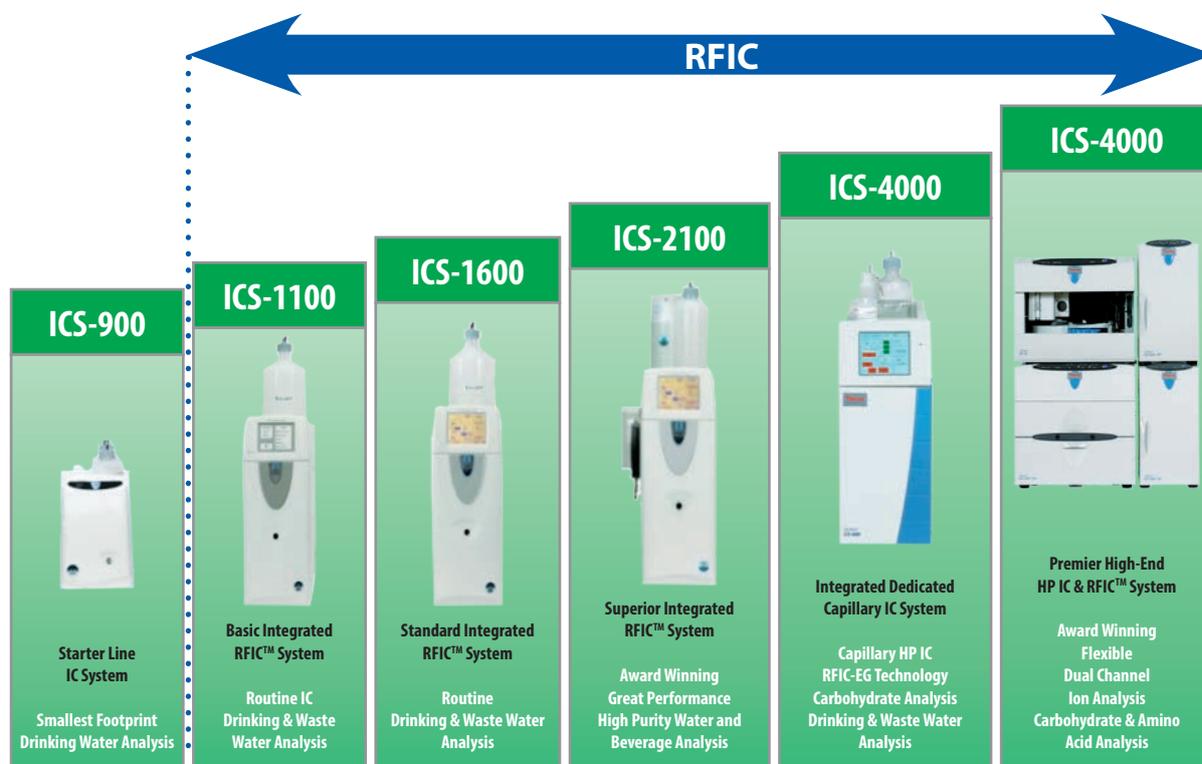
ION CHROMATOGRAPHY

Thermo Scientific™ Dionex™ ion chromatography (IC) systems have led the industry for over 30 years, with instruments and solutions that represent state-of-the-art technological advancements and patented technologies, including:

- High-Pressure IC
- RFIC™ System technology (Reagent Free IC)
- Eluent generation
- Electrolytic suppression
- Capillary IC
- Continuously regenerated trap columns
- Three-dimensional electrochemical detection
- Award-winning systems
- Dual capabilities to improve workflow, signal enhancement, and complex applications

A Wide Range of Solutions

Thermo Fisher Scientific develops, manufactures, sells, and services IC systems to separate, isolate, and identify the components of chemical mixtures. Dionex Products IC systems and separation chemistries are used extensively in the environmental, food and beverage, pharmaceutical, life science, biotechnology, chemical, petrochemical, power generation, and electronics industries. Applications include anions and cations, carbohydrates, organic acids, amino acids, proteins, peptides, oligonucleotides, oxyhalides, carboxylates, polysaccharides, polyphosphates, fatty acids, metals, phenols, surfactants, ethanolamines, and more.

Systems

Depending on your needs, a wide range of choices are available, from the Thermo Scientific Dionex ICS-900 to the Dionex ICS-5000+ HPIC™ system. Whatever the size of your workload, whether your analytical task is simple or challenging, we have a solution to match your needs and budget. Dionex IC systems also come with superior customer support and service worldwide.

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ATOMIC ABSORPTION SPECTROSCOPY

The method of atomic absorption spectrometry (AAS) allows determining a wide range of elements: metal ions (alkali and alkaline earth), heavy metals, toxic elements in the water and in the soil.

240 AA are atomic absorption spectrometers with easy-to-use software at a mid-range price.

240FS AA fast sequential atomic absorption spectrometer brings the productivity and analysis speed of ICP-OES to atomic absorption.

240Z AA (4-lamp) offers the highest sensitivity, performance, simplicity of operation, and industry-leading software in atomic spectroscopy.

280FS AA flame instrument combines Agilent's Fast Sequential design with 8-lamp capability for greater performance and reduced running costs.

280Z AA (8-lamp) offers the highest sensitivity, performance, simplicity of operation and industry-leading software in atomic spectroscopy.

SpectrAA 55 series is perfect for mobile labs in remote location or any laboratory needing an entry-level Atomic Absorption system.

AA Duo is the world's only atomic absorption spectrometer with simultaneous operation of flame and graphite furnace.

Methods for water:

- GOST 51309-99 Drinking water. Determination of elements by atomic spectrometry.
- GOST 51212-98 Drinking water. Methods for determination of total mercury by flameless atomic absorption spectrometry.
- АРНА 3111 Metals by flame atomic absorption spectrometry.

Methods for soils:

М-МВИ-80-2008 Methods for measuring the mass fraction of elements in soil samples, soil and sediments by atomic emission and atomic absorption spectrometry.

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UV-VIS-SPECTROSCOPY

Cary 60 UV-Vis Spectrophotometer

The Cary 60 UV-Vis Spectrophotometer is efficient, accurate and flexible, and is designed to meet your immediate and future challenges.

Measure precise samples with ease – the highly focused beam image of the Agilent Cary 60 is perfect for measuring small volumes accurately and reproducibly. Preserve your samples by using < 4 µL instead of mLs

Exceptionally fast data collection – with a scan rate of up to 24,000 nm/min, you can scan the entire wavelength range (190-1100 nm) in under 3 seconds

Performance enhancing accessories – assures you can handle the widest variety of sample sizes and types with liquid sampling accessories including fiber optic probes and couplers, Peltier and water thermostatted single and multicell holders, temperature probes, microvolume sampling cells and rapid mix accessory for stopped flow kinetics. For solid samples, choose from solid sample holder, fiber optic reflectance probe and coupler and fixed angle specular reflectance accessories (SRA)

Software designed for real samples – the Agilent Cary WinUV software can be tailored to suit your analytical requirements from QA/QC to research

Specification	
Monochromator	Czerny-Turner
Detector	Twosiliconphotodiode
Diffractiongrating	Holographic, 27.5x 35mm, 1200lines / mm, blazeangle – 8.6 °at 240 nm
Spectralrange	190-1100nm
Spectralslitwidth	Fixed, 1.5 nm
Wavelengthreproducibility	± 0.1nm
Wavelengthaccuracy	± 0.5nm
Max. scanspeed	24,000nm/min

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DR Series Spectrophotometers

Upscale spectrophotometers turnkey solutions for both portable (DR-2800, DR-3900) and laboratory (DR-6000) analysis. Designed specifically for water analysis. Allows to work with process waters and with drinking or waste water.

Application areas:

- Municipal and industrial wastewater;
- Drinking water and surface water;
- Water treatment;
- Production and quality control.

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DR/6000 - professional laboratory UV / VIS spectrometer

The DR6000™ is advanced laboratory spectrophotometer. It offers high-speed wavelength scanning across the UV and Visible Spectrum, and comes with over 250 pre-programmed methods including the most common testing methods used today. With optional accessories allowing for high-volume testing via a carousel sample changer, and increased accuracy with a sample delivery system that eliminates optical difference errors, this instrument ensures you are ready to handle your wide-ranging water testing needs.

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Specification	
WavelengthRange	190...1100 nm
WavelengthAccuracy	±1 nm (200...900 nm)
WavelengthReproducibility	< 0.1 nm
WavelengthResolution	0.1 nm
Scan Speed	900 nm/min. (in 1 nm step)
Spectralbandwidth	2 nm
Photometric accuracy	±0.005 Abs (0.0...0.5 Abs); ±1 % (0.5...2.0 Abs)
Photometric linearity	< 0.5 % (<2 Abs); < 1 % (>2 Abs)
Light scattering	< 0.05 %

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HIGH PERFORMANCE LIQUID CHROMATOGRAPHY

High performance liquid chromatography (HPLC) is one of the most effective methods for separating complex mixtures of compounds, widely used in analytical chemistry as well as in chemical technology.

UltiMate 3000 series intelligent LC (LCi) HPLC systems are one of the world's best developments for high-performance liquid chromatography for simple and reliable operations. A wide selection of modules allows you to optimize the system for your specific purposes: from ultra-high performance rapid separation liquid chromatography (RSLC) to highly sensitive nano LC / MS.

Whichever configuration you choose, you get a complete system with optimized fluid paths and high-performance software Chromeleon® for the management, control and data processing.

UltiMate 3000 Intelligent LC System won the gold medal of Instrument Business Outlook Journal on the Pittsburgh Conference as the best product of the year.

Key characteristics and features of the system

- Rapid separation liquid chromatography (RSLC)
- Rapid separation columns for RSLC
- Excellent quality of chromatographic separation
- Unique dual systems for high performance and complex procedures
- Reducing of analysis time, speeding up the process of validation
- Automated many dimensional separation
- System with switching of automatic methods
- Ease of use

Variety of systems

- Rapid separation liquid chromatography (RSLC)
- UltiMate3000 HPLC system in the titanium version, designed specifically for the analysis of biomolecules.
- Classic system
- Quaternary analytical system
- The binary analytical system
- Isocratic analytical system
- Basic manual and automatic systems
- Micro/nano system
- Quaternary micro system
- The binary micro system
- Quaternary nano, capillary and micro systems

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ORGANIC COMPOUNDS ANALYSERS

DRB200 Digital Reactor for COD of water

DRB200 Digital Reactor for analysis of COD and cuvette tests. COD ranges 3–150 mg/L and 20–1500 mg/L COD are USEPA approved (5220 D) for wastewater analyses.

Decomposition can be accomplished using the potassium permanganate method, as well as the dichromate method. After decomposition of the sample, the measurement is made using a spectrophotometer DR series (3900, 6000) or other UV/Vis-spectrophotometers.

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BOD analyser OxiTop® IS 6, IS 12 with OxiTop® Box incubator APHA 5210 Biochemical Oxygen Demand (BOD)/ D Respirometric Method

Complete packages for 6 or 12 simultaneous measurements. Measurement using OxiTop® is based on pressure measurement in a closed system: microorganisms in the sample consume the oxygen and form CO₂; the CO₂ is absorbed by NaOH, creating a vacuum that can be measured as a mg/l BOD value. The sample volume used regulates the amount of oxygen available for a complete BOD. Measurement ranges of up to 4,000 mg/l can be measured using different volumes. The OxiTop® heads (green and yellow for inflow/outflow differentiation) have an AutoTemp function: if the sample temperature is too cold, the start of measurement is automatically delayed by at least 1 hour until a constant temperature has been reached. Apart from the automatic storage of 5 measured values (1 value per day), further measured values can be read at all times during or after the period of 5 days, which permits the tracking of check values or measurements over longer periods.

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ANALYSERS OF OIL PRODUCTS IN WATER AND SOILS

OCMA-350 Oil Content Analyzer

Compact and easy to use, the OCMA-350 oil content analyzer provides quick, accurate measurements in the lab or in the field. The instrument measures the amount of oil in water, soil and machined parts for a variety of applications. Some common applications include waste water effluent, bilge water discharge, oil/water separators, contaminated soil and oil on textiles and metal parts. The OCMA-350 uses an excellent extraction solvent called S-316.

Features

- The instrument measurement takes 30-60 seconds using a solvent for the extraction and NDIR method.
- The use of a quartz sample cell keeps the instrument clean from possible spills.
- Oil is extracted using the extremely stable S-316 solvent extractant.
- Users can display analysis results in three different units-mg/l, mg/kg, and Abs.
- The instrument has an easy-to-read backlit liquid crystal display.
- RS-232C interface is standard equipment.

Recommended for application in the environmental monitoring, analysis of water quality, wastewater analysis, in the marine transportation, refining, monitoring on tank farms, etc.

ASTM D7066 - 04(2011) Standard Test Method for dimer/trimer of chlorotrifluoroethylene (S-316) Recoverable Oil and Grease and Nonpolar Material by Infrared Determination

GOST R 54039-2010 Soil quality. Express-IR spectroscopy to determine the number and identification of soil contamination with oil products

PND F 16.1:2.2.22-98 method for measuring the mass fraction of oil in mineral, organogenic, organo-mineral soils and sediments by infrared spectrometry

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TOTAL ORGANIC CARBON (TOC) OF WATER AND SOIL

Vario TOC cube

Elemental analyzer **vario TOC cube** designed for the simultaneous determination of total organic carbon and other forms of carbon (total carbon, total inorganic carbon, dissolved carbon, volatile and non-volatile organic carbon), both in solid and in liquid samples or suspensions.

Specifications

Measuring method

Unlike other methods of decomposition, the method of high-temperature combustion in oxygen allows more complete decomposition of the organic compounds present in the sample and get more accurate results. High-temperature furnace provides heating up to 1200 °C. Lifetime warranty of the oven - 10 years!

Large detection range

0 to 60,000 mg/L with a detection limit of 6 ppb. Variable sample volume between 0.05 and 2 mL allows also multipoint calibration from a single standard solution with auto-dilution-effect.

Modular concept for liquids and solids analysis

Large choice of automatic sample feeders for liquids and solids, alternatively cost-efficient solutions with manual sample change. With possibility to upgrade at any time.

Real matrix separation

The new reactor concept enables real matrix separation and offers maximum protection of the combustion tube. Excellent properties for analyses of saline solutions (e.g. brine) or other complicated sample matrices.

Software with high performance and maximum flexibility

Windows® XP professional or Vista business based control and evaluation software with calibration wizard, system suitability test, statistics and advanced math operations. Integrated maintenance software and comprehensive help menu.

Supports validation

Complete qualification documentation with Standard Operation Procedures (SOPs) and Certificates for IQ, OQ and PQ according to FDA regulations

Application

GOST R 52991-2008 Water. Methods for determination of total and dissolved organic carbon
GOST 17.1.5.03-81 Conservancy. Hydrosphere. Total organic carbon analyzers in natural waters

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EQUIPMENT FOR AIR QUALITY MONITORING

For air quality monitoring both atmospheric and working area can be used miscellaneous equipment - gas analyzers both stationary and portable, aspirators, indicator tubes, sampling devices and filters.

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STATIONARY GAS ANALYSERS

T-Series analysers, manufactured by Teledyne-API, can be at monitoring of air quality, continuous emissions monitoring and control of indoor air with very clean atmosphere.

Common features of all T-Series gas analyzers:

- Large, vivid, and durable color graphics display with touch screen interface
- Front panel USB connections for peripheral devices and firmware upgrades
- Internal data collection system for storing data and diagnostic messages
- Ethernet, RS-232, and (optional) USB - COM ports
- Common line of spare parts suitable for many gas analyzers

Physical Specifications:

Dimensions (D x W x H), mm: 597 x 432 x 178

Weight, kg: from 18 to 30 (depending on the specification and configuration of the device)

Operating Temperature Range: +5°C ... +40°C

To calibrate the analysers can be used as built-in calibration sources on the basis of microflow sources - (optional), as using external calibration sources.

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STATIONARY GAS ANALYSERS

Gas	Model	Ranges (min / max)
Measuring principle: Ultraviolet Fluorescence		
SO ₂ Sulfur Dioxide	T100U	0...5ppb / 0...20ppm
	T100	0...50ppb / 0...20ppm
	T100H	0...10ppm / 0...5000ppm
H ₂ S Hydrogen Sulfur	T101	0...50ppb / 0...20ppm
TRS Total Reduced Sulfur	T102	0...50ppb / 0...10ppm
TS Total Sulfur	T108U	0...5ppb / 0...20ppm
	T108	0...50ppb / 0...20ppm
Measuring principle: Chemiluminescence		
NO, NO ₂ , NO _x Nitrogen oxide (II) Nitrogen dioxide (IV)	T200U	0...5ppb / 0...2ppm
	T200PU (uses photolytic converter)	0...5ppb / 0...2ppm
	T200	0...50ppb / 0...20ppm
	T200M	0...1ppm / 0...200ppm
	T200H	0...5ppm / 0...5000ppm
NO _y Oxides of Nitrogen	T200U / NO _y	0...5ppb / 0...2ppm
NH ₃ Ammonia	T201	0...50ppb / 0...2ppm
O ₃ Ozone	T265	0...100ppb / 0...2ppm
Measuring principle: Infrared Gas Filter Correlation		
CO Carbon Monoxide	T300U	0...100ppb / 0...100ppm
	T300	0...1ppm / 0...1000ppm
	T300M	0...5ppm / 0...5000ppm
CO ₂ Carbon Dioxide	T360U	0...100ppb / 0...100ppm
	T360	0...2ppm / 0...2000ppm
	T360M	0...4ppm / 0...4000ppm
N ₂ O Nitrous oxide	T320U	0...200ppb / 0...200ppm
	T320	0...1ppm / 0...1000ppm
Measuring principle: Ultraviolet Absorption		
O ₃ Ozone	T400	0...100ppb / 0...10ppm
Measuring principle: CAPS (cavity attenuated phase shift)		
NO ₂ Nitrogen dioxide (IV)	T500	0...5ppb / 0...1ppm

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EQUIPMENT FOR CALIBRATION OF GAS ANALYSERS

To ensure uninterrupted operation with the specified accuracy characteristics analyzers must be calibrated. For the calibration can be used as calibration gas mixtures with specified concentrations, as sources of the microflow.

Gas	Model	Ranges (min / max)
Dynamic Dilution Calibrator	T700U	0...100%
	T700	0...100%
	T700H	0...100%
Air Zero Generator	701	0...20 SLPM, 30 psi
	701H	0...30 SLPM, 35 psi
Portable Air Zero Generator	751H	0...10 SLPM, 30 psi
	751	0...5 SLPM, 30 psi
Ozone (O3) Calibrator	T703U	Ozone concentration: 3ppb...5ppm,
	T703	Ozone concentration: 50ppb...5ppm

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Dynamic Dilution Calibrator T700

Dynamic Dilution Calibrator model T700 allows the dilution of the calibration gas mixture ranging from 0 to 100% of the concentration. As a diluent gas used "zero air" from a cylinder or zero air generator. Calibrator can simultaneously connect up to 4 cylinders with single or multi-component gas mixture. Calibrator-diluent may be equipped with UV ozone generator for calibration of ozone analyzers and test drive NO to NO₂. For more precise control of generating ozone calibrator-diluent may be fitted with a photometer.

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Zero Air Generator 701

Zero Air Generator produces dried and free of impurities air - «zero air». Depending on the application, the generator can be equipped with filters to remove carbon oxides and hydrocarbons

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Portable zero air generator T751

Makes it possible to produce «zero air» as the model 701, but in small quantities.

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PORTABLE GAS ANALYSERS

Portable flue gas analyser Testo 350

Testo 350 is a flexible portable flue gas measuring system wide with range of applications for combustion systems. The equipment can be used for:

- For all types of industrial boilers settings
- Long time measuring of exhaust gas concentrations
- Operational control of industrial burners
- Monitoring and adjustment of stationary diesel generators
- Measurement of pressure and velocity in ducts and flues

Analyser TESTO 350 is designed to control the flue gases and can measure O_2 (factory setting) and the concentration of the following gases: CO , CO_{low} , NO , NO_{low} , NO_2 , SO_2 , H_2S , HF , CO_2 (separate IR sensor, or calculation), the calculation of NO_x , and flue gas flow rate (using the Pitot) performance factor, traction / pressure built option differential pressure measurement, a trigger input, USB-output, probe connector for temperature measuring (type K NiCr-Ni and type S Pt10Rh -Pt).

TESTO 350 control unit is designed to control the measuring system and display of measurement data. The module can be removed from the analyser and upgraded with lithium-ion battery. All settings are made using the arrow keys. The measurement results are displayed on the color graphics display. Built-in memory data can be transferred from the analyser to the control unit. It is possible to convenience control several analyser unit by single control unit.

- Remote control of analyser unit and measurement data transfer from chimney flue. It's very convenient for industrial burner measurement.
- Measurement data transfer from analyser unit to control unit allow to remote control unit from analyser and retrieve obtained data from analysers. The analyser unit keeps operating at that time.

To protect the display of control unit Testo 350 from being damaged by long-term measurements or transportation, the control unit can be installed in the analyzer unit in face down position.

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Portable flue gas analyser Testo 340

Compact shape and available equipment make analyser testo 340 ideal for operation, service and maintenance, as well as control measures of industrial boilers, gas turbines, stationary industrial engines, thermal processes.

Advantages of analyser Testo 340:

- Maximum quantity by sensors: 4.
- The standard unit is equipped with O_2 sensor. 3 additional sensors be can configure by your own choice: CO (H_2 -compensated), SO_{low} (H_2 -compensation), NO, NO_{low} , NO_2 and SO_2 .
- Standard built-in measuring system of the differential pressure and velocity.
- CO (H_2 -compensated) and NO sensor with replaceable filters. This filter prevents the gas passage through the sensor. If the filter is worn out, the user can replace it by himcell.
- Optimum shockproof protection.

Testo 340 – calculated parameters

- dew point
- heat loss +performance
- CO – undiluted
- NO_x
- flow rate (volume flow, mass flow)
- λ (lamda)
- gas consumption
- battery capacity (%)
- sensor diagnostics

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GANK-4 gas analyser

GANK-4 analyser A, R, AR modification designed for automatic continuous monitoring of concentrations of harmful substances in the air (A), in the working area (R), industrial emissions and processes for environmental protection, occupational safety and process optimization

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GANK-4 CONTROLS 150 HARMFUL SUBSTANCES FOR CHOICE IN THE MEASUREN POINT WITHOUT SAMPLE PREPARATION

Specification:	
Range of the GANK-4(A)	from 0,5 MPC d.a.* to 0,5 MPC w.z.**
Range of the GANK-4(R)	from 0,5 MPC w.z.** to 20 MPC w.z.**
Range of the GANK-4 (AR)	from 0,5 MPC d.a.* to 20 MPC w.z.**
Measurement time	10-30s.
Measurement error, no more	20%
Dimensions, mm	250x200x150
Weight ,kg	Not more 3,5
Power consumption, W	Not more 8
Ambient temperature	From - 50°С to + 50°С
with thermostat TP-1	From - 50°С
Internal battery	12V
Mains power	220V, 50Hz

Various components are measured exchangeable cassette or built-in sensors. Each cassette contains an electronic memory that recorded measured substance and measurement parameters.

* MPC d.a. - Average daily maximum permissible impurity concentration in the atmosphere, in mg/m³ (for substances which are produced by the sensors).

** MPC w.z. – The maximum permissible concentration of harmful substances in the air of the working area, in mg/m³.

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HIGH VOLUME SAMPLERS

High Volume Samplers DHA-80.

High Volume Samplers DHA-80 are fully automatic systems to sample dust and aerosol particles for later assessment and analysis (gravimetric and analytical determination).

The sampler operation range in standard execution is 100 to 1.000 liters per minute (6 to 60 m³/h).

Sampling is done on the filters installed in the sampler. 15 liters can be stored in the sampler.

Operating principle: switching on, sample tacking, switching off, pausing and cycle repeat.

Sample taking is from 0 to 60 minute including pause.

Dimensions (L x W x H): 600mm x 250mm x 1385mm (excluding sampling probe).

Weight: depending on the specification and configuration up to 60kg.

Operating temperature range: from +5 ° C to +40 ° C and humidity from 10 to 90% or -20 ° C to +40 ° C and humidity of 10 - 95% with built-in heating system.

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WIND SPEED AND WIND DIRECTION SENSORS

Combined sensor model 034B for wind speed and wind direction measurement.

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Model 50.5 ultrasonic sensor wind speed and wind direction.

Depending on the operating conditions the model is supplied with standard and extended temperature ranges.

Features:

- Without moving parts
- Analog and digital outputs
- Can be supplied with heating circuits that greatly extend the temperature range during of operation

Specifications:

- wind speed:

- Range: 0 ... 50 m/s
- Accuracy 0,1 m/s
- Resolution 0,1 m/s

- wind direction

- Range: mechanical 0 – 360°
- Accuracy $\pm 3^\circ$
- Resolution 1°

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BRIEF OPERATION DESCRIPTION OF AQMS (AIR QUALITY MONITORING STATION)

Air quality monitoring station (AQMS) represents a stationary enclosure with dimensions 4x2,5x2,8m, which may vary depending on the quantity of primary and secondary equipment, with telescopic mast for meteorological equipment and area for maintenance sampler system.

Depending on customer requirements enclosure can be installed either on the ground or concrete base, and the metal pedestal.

AQMS provides continuous control of ambient air for determining air pollutants.

Basic version of AQMS configurations defines following air pollutants: mercaptans, $\text{NO}/\text{NO}_2/\text{NO}_x$, NH_3 , O_3 , CO_2 , CO , SO_2 and H_2S .

List of air pollutants may be vary and agreed with the Customer.

Measuring of air pollutants in ambient air are carried out continuously using gas analyzers.

Air is injected by the PTFE tube with the help of the pump to sampler manifold. From the manifold the air enters to the analyzers by the PTFE tubes.

Excess air is extracted out.

In addition to measuring of the concentration of air pollutants by appropriate sensors it is monitored such meteorological parameters as wind speed and wind direction, barometric pressure, humidity and air temperature.

Every 2 seconds data is taken from analysers and sensors and using input modules converted and transmitted to the PC.

Measured data on both concentration of air pollutants, and meteorological parameters enters to analog input modules.

IO module of discrete signals controls the analysers status, as well as an operation of enclosure heater, sampler pump and door status.

A software is installed in PC which performs remote monitoring and controls of devices and actuating mechanisms as well as it displays on the monitor the current value of all devices and indicates the maximum permissible con-

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centration (MPC) of substance in air, the calculation of the average values within the time interval (agreed with the Customer). All data is stored in the database.

Life-support system includes two heater and air conditioner to maintain the optimal temperature range for normal operation range of devices.

Indoor temperature is controlled via temperature sensor installed on rack with devices.

Also fire and secure alarm system is provided in the AQMS. Upon the customer request the information from fire and secure alarm system can be transmitted to the customer's security service, i.e. the company responsible for protection of Customer's objects.

Power supply of AQMS provided from AC line 220V or 380V via UPS. All equipment installed in the AQMS has an operating voltage of 220V, 50Hz.

In case of power failure the battery life of the uninterruptible power supply of the AQMS is determined by the Customer. Backup power is performed either from additional power lines or from a backup generator, or may not be available. If you intend to install several such AQMS at the facility for simultaneous and continuous air monitoring at several points in the necessary area.

AQMS will be connected with central monitoring station for transmitting data of average value. Received from stations data results can be transmitted both by radio, cable lines, and via satellite terminal, depending on customer preference.

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